


# T-Bot® II Class Pack

## User Guide

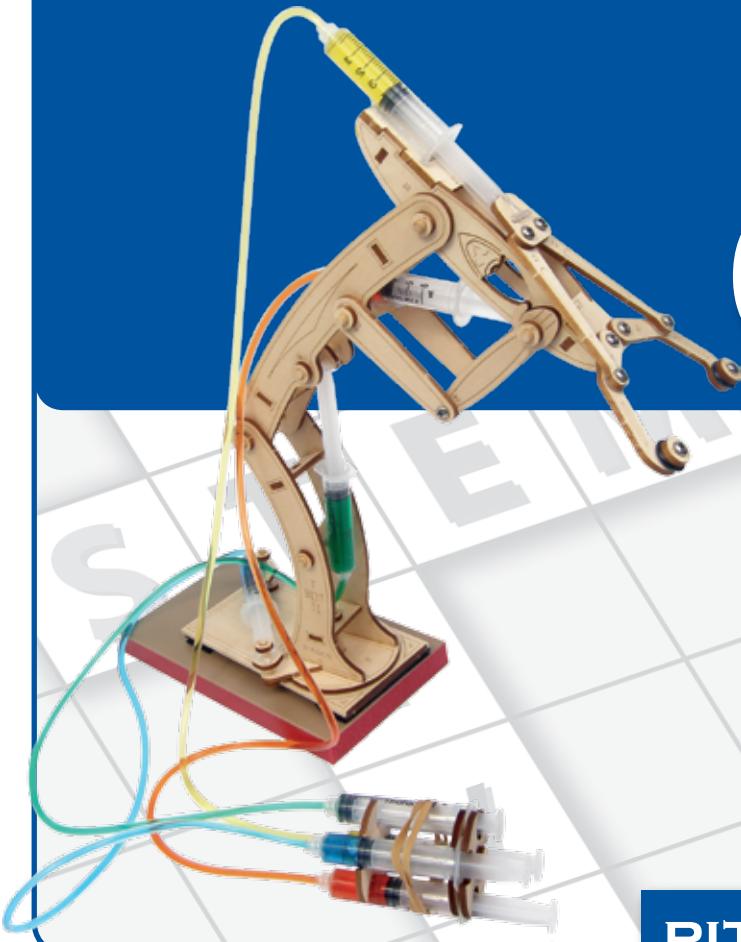

**PITSCO**  
EDUCATION

The T-Bot® II Class Pack is bulk-packed to save you money. As such, you need to divide the class pack into 10 individual T-Bot II kits. To do this, follow the instructions below. If anything is missing from your class pack, please contact Customer Service at 800-358-4983.

1. Remove all items from the box.
2. From the plastic bags, remove the basswood sheets labeled with a letter "A." Spread these out on a large, clear surface such as a table or counter.

**Note:** If any laser-cut parts fall out of a sheet, make sure to keep these parts in the correct pile so no kit is missing parts.

3. From the plastic bags, remove the basswood sheets labeled with a letter "B." Place one with each of the Sheet 1 basswood pieces.
4. From the plastic bags, remove the smaller basswood sheets labeled with a letter "C." Place one with each of the Sheet 1 and 2 basswood pieces.
5. Continue this process with the sheets labeled with the letters "D" and "E."
6. Place one Precision Straw and one resealable bag on each pile.
7. Add one rubber band, two 1/2" screws, and eight 3/8" screws to each pile.

*(continued on back)*

8. Place two O-rings and two nylon spacers on each pile.
9. To each pile, add the following number of dowels:
  - 1 – 1-1/8" dowel
  - 1 – 1-1/2" dowel
  - 4 – 1-3/4" dowels
  - 3 – 2" dowels
  - 1 – 2-1/2" dowel
10. Place eight syringes in each pile.
11. Find the length of clear tubing. Using scissors, cut the tubing into eight-foot lengths until you have an eight-foot length for each pile. You will have approximately 20 feet left over.
12. There will be 5 one-foot lengths that are the two sides of the hook-and-loop fastener. Cut each of these lengths in half to make 10 pieces (of each side of the fastener). Add one piece to each pile.
13. Finally, add one T-Bot II User Guide to each pile.
14. Place each pile in a bag, box, or some other container to keep the parts together until you use them in class.



P.O. Box 1708 • Pittsburg, KS 66762  
shop.pitsco.com • Toll-Free Orders 800-835-0686  
For Customer Service,  
call 800-358-4983 or e-mail [service@pitsco.com](mailto:service@pitsco.com)